

TML\MSH Microbiology Department Policy & Procedure Manual	Policy # MIMYC\09\12\v01	Page 1 of 1
Section: Mycology Bench Manual	Subject Title: Media / Reagents Sabouraud Gentamicin Agar Slopes	
Issued by: LABORATORY MANAGER	Original Date: March 22, 2001	
Approved by: Laboratory Director	Revision Date:	

MEDIA / REAGENTS

12. SABOURAUD GENTAMICIN (50 mg/L) AGAR SLOPES

Purpose

To isolate fungi especially yeast from contaminated specimens.

Sabouraud Agar (Difco-modified-0747-01)	50 g.
Dist. H ₂ O	1000 ml.

Mix well, bring to a boil to dissolve. Cool to 50⁰C (check pH with ATC - pH ± 7.0). Add 1.25 ml Gentamicin (1 vial = 80 mgm/2 ml.). Mix well. Dispense 10-ml amounts into pre-sterilised UGB bottles. Autoclave 118⁰C/10 minutes. Cool in a slanted position. Tighten the caps. Label the bottles "SAB G". Store at 4⁰C.